




Switching Between VCOM and USB Keyboard for CODE CR1500 Scanners

The instructions below will allow you to switch between USB Keyboard and VCOM settings. If you haven't done so already, please make sure that you have properly integrated your scanner first. More information may be found on the RapID Lab website:

<http://rapidlab.com/scanner-and-software-support/>

USB Keyboard Settings:	VCOM Settings:
<p>1. Scan the following barcode to remove any previous programming from the reader</p>  <p>CC004361_02</p> <p>2. Scan the following barcode to program the reader for custom settings requiring a suffix<ENTER></p>  <p>CC004482_03</p>	<p>1. Scan the following barcode for enabling VCOM and disabling data formatting</p>  <p>CC004854</p>

Additional optional codes available on the following page.

Optional Codes:

1. Scan the following barcode to disable suffix <ENTER>



M20208_01

2. Continuous Scan On



M20224_01

3. 1 Second Duplicate Scan Delay



M20230_01

4. 3 Second Duplicate Scan Delay



M20232_01

1. This code disables the suffix <ENTER>.

2. This code allows for continuous scanning on your CR1500.

3. This code creates a 1 second delay between scans.

4. This code creates a 3 second delay between scans.

5. Disable Duplicate Scan Delay



M20229_01

5. This code disables duplicate scanning.