

## Setting up CODE CR1500 Scanner for Rapid Tags as a USB keyboard

The instructions below will program the CODE CR1500 Scanner to properly decode the Rapid Tag Data matrix symbols. More information may be found on the Rapid Lab website:

<http://rapidlab.com/scanner-and-software-support/>

1. First ensure that there is no plastic packaging/protective film attached to the scanner lens.
2. Next, plug in the scanner to your computer, laptop, or tablet and open a simple word document or notepad for testing.
3. Scan the below barcodes in the given order. When scanning the configuration codes, briefly hover the light on top of the barcodes and quickly remove to ensure that each code is scanned only once.

### Required Codes:

1. Remove previous programming (you should hear 3-4 beeps)



CC004361\_02

2. Custom Scanner Settings



CC004482

4. Next, test your new scanner configurations by scanning a Rapid Tag which should populate with a 6 digit alphanumeric ID and a line break after.