



Rapid Lab Recommended Sterilization and Disinfecting Procedures

Please note that RapID Lab products do not ship sterilized. We highly recommend sterilizing and/or disinfecting your equipment prior to use.

Please refer to the steps below for the best results:

1. Remove your RapID Lab products from the packaging and place the items in an autoclave bag.
2. Run at a cycle of 121deg Celsius (249deg Fahrenheit) for 25 minutes and let it cool completely before removing for use.
3. Before applying the tags, give your applicator and tags a quick rinse with 70% isopropyl alcohol to wash off any potential dust or debris then let dry.